


Quick Start Guide

Flexible Wavelength Selector - Auto



1

Connect the power supply of FWS Auto to a power outlet

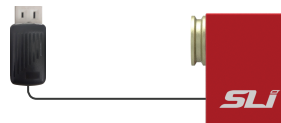

2

The ►►► arrow on the image indicates the directionality of the propagating light.




3

Download and install the FWS Auto software from www.spectrolightinc.com. Connect the FWS Auto to a computer using the provided USB cable.



4

The center wavelength and the FWHM can be selected using the SET of the provided software.

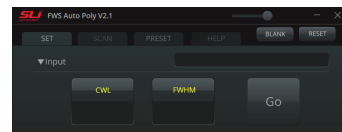

5

The SCAN tab provides options to scan a range of wavelength with user selected FWHM, step size and time delay.



✳ For more details, please refer to our website www.spectrolightinc.com