

Quick Start Guide

Flexible Wavelength Selector - CenterLine



1

▶▶▶ This arrow indicates the directionality of the light source.



2

As seen on the image at the right, there are two knobs. The one on the left (1) can be used to select your desired wavelength while the one on the right (2) is used to precisely compensate the directionality offset of the light.



3

On the knobs you will be able to see markings L and S which refers to long and short wavelength, respectively.



4

By referring to the Wavelength Guide Table in our website, it is possible to select the center wavelength. For more accurate determination of the center wavelength and the bandwidth, please use a spectrometer.



※ For more details, please refer to our website www.spectrolightinc.com